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Control panel

The control panel is fully digitised and contains all the operating controls and the panel printer.

Information regarding all the cycle parameters, time, temperature and pressure are displayed on specific displays of the control panel to allow clear reading at a glance.

There are no manometer or analogue controls or systems.

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Preheating system

The preheating system has been designed to be quick and to guarantee rapid, controlled cycles.

When the equipment is not in use, a special system allows the chamber to be kept in thermal stand by mode, so guaranteeing reduced preheating times.

After an hour of non-use the equipment enters power saving mode and switches itself off.

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Errors and alarms

The autoclave has an integrated control system that is able to detect technical anomalies and shifting of cycle parameters, which it indicates with ALARM messages.

The check detects ERROR situations caused by incorrect use or incorrect operations.

In this way, the operator can have a continuous diagnosis of the equipment and be certain that it is in working order.

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Sterilization chamber

The sterilization chamber is made of AISI 304 stainless steel.

It has a diameter of 245 mm, is 500 mm deep and can be provided with a basket.

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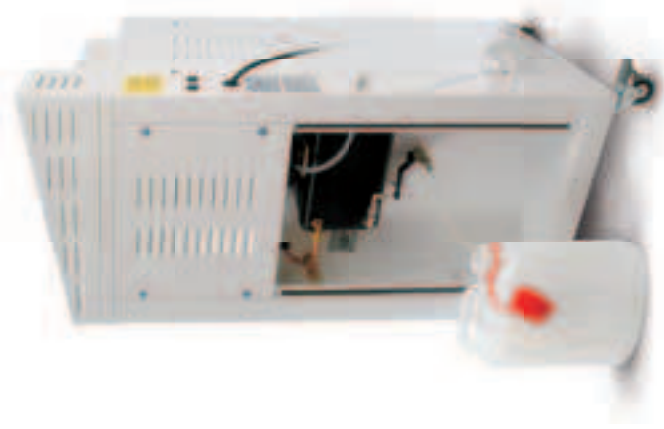
LABCLAVE has a wide range of programs that are also suitable for sterilizing liquids.

There are 4 automatic programs set with different decompression systems; rapid or normal and 1 free cycle that can be programmed according to the operator's requirements.

LABCLAVE also has a system called NIGHT CYCLE that is activated automatically on all the programs 60 minutes after the end of the cycle and automatically switches off the equipment.

To assess the response of the cycle, the printer document can be assessed or any control button can be pressed to make the machine automatically switch on the displays that will have held any alarms or messages in the memory.

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Tank

The autoclave carries out **automatic draining of recondensed steam** into a special container with overflow LED indicator.

LABCLAVE is also provided with a tap to enable draining of residual water or water present inside the sterilization chamber.

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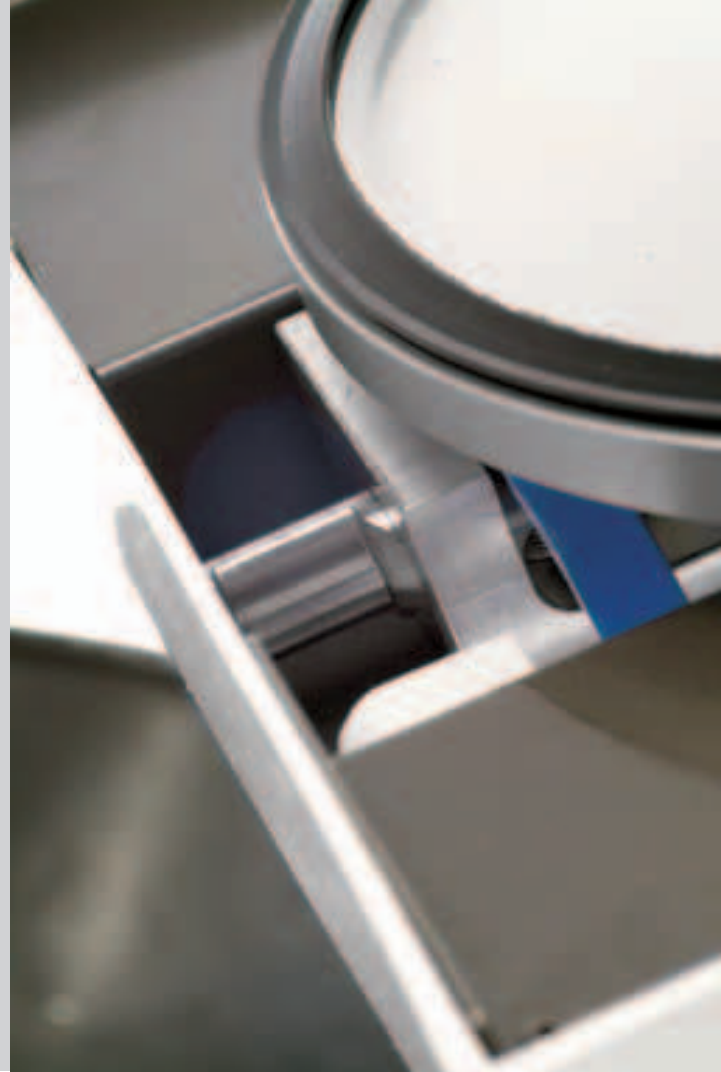
The autoclave is manufactured according to European standards and rigorously applies all the safety systems.

LABCLAVE is equipped with:

- certified safety valve for overpressure management
- chamber insulation with certified ceramic material

It also has a safety system on the door that prevents the possibility of it being opened in the presence of internal pressure.

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Close-up view of the front of an Labclave

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Tecno-Gaz uses all the standards required by safety legislation in force. But in addition to that we have decided to comply with these regulations adding also other technical solutions for our customers.